


# Tk Novafix™ HPLC Cartridge System

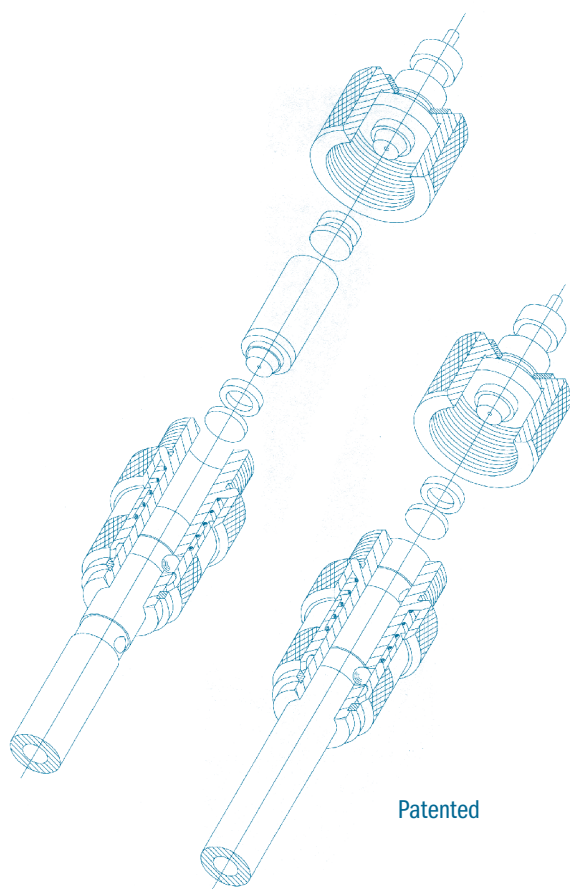

Teknokroma has designed and developed the original, patented Novafix™ Cartridge System for HPLC, which is the result of more than 14 years of experience researching and manufacturing Teknokroma HPLC columns.

## Novafix™ Quality

Teknokroma has always achieved its best in offering top-quality products and services. This quality-excellence philosophy has helped Teknokroma achieve HPLC market leadership wherever its products are marketed. Our research scientists have utilized the same quality-excellence philosophy to meticulously design the new Novafix™ HPLC Cartridge System. Novafix™ Cartridges are made from chosen materials and select bonded packings that guarantee the greatest column efficiency, peak symmetry and reproducibility.

In addition, Teknokroma's proprietary Novabond™ packing procedures are the result of years of exhaustive research and detailed manufacturing of HPLC columns. Novabond packing procedures provide you with the best column efficiency, peak symmetry and column lifetime available on the market.

## Novafix™ Easy Handling

The mechanism for rapid connection designed in the Novafix™ Cartridge System does not require you to use any tools for its assembly or dismantling. This design feature makes Novafix™ columns simple and easy to handle.

## Novafix™ Functional Design

Without requiring any additional accessories, the Novafix Cartridge System permits the insertion a 1cm-long precolumn at the head of the analytical cartridge. This is achieved without introducing any dead volume, thereby maximizing column efficiency and peak symmetry.


## Novafix™ Efficiency

The Novafix™ HPLC Cartridge supplies the highest theoretical efficiency. These typically high efficiencies are achieved due to the zero dead-volume connections and proprietary Novabond™ packing procedures.

Particle size $\mu\text{m}$	Typical Efficiencies N/m
3	120-150.000
5	80-110.000
10	35-65.000

### Typical Efficiencies

Efficiency media: 85300  
St: 5298

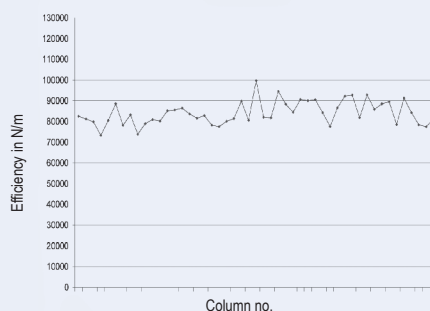


ODS type columns.  
Efficiencies calculated with the peak of Anthracene using ACN/H<sub>2</sub>O 65:35 as eluant.

## Novafix™ Stability

The design of the Novafix™ HPLC Cartridge System ensures not only maximum efficiencies, but also long useful lifetimes. Novafix™ Cartridges provide maximum stability for packing materials in the precolumn and analytical column cartridges, no matter how frequently the precolumn is exchanged.

### Stability

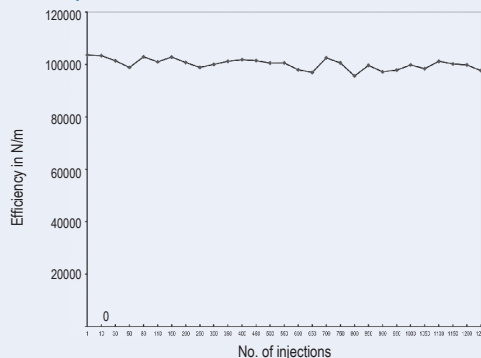


# Novafix™ HPLC Cartridge System Tk



## Novafix™ Reproducibility

In the Novafix™ cartridges manufacture, we only use the top valued packings of the market for guaranteeing the maximum reproducibility. That way the values of resolution, selectivity, efficacy and stability will repeat column by column and year by year, without requiring special modifications in the chromatographic method.

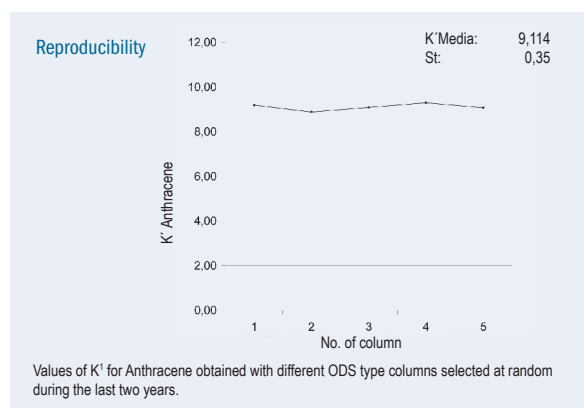

## Novafix™ Versatility

Teknokroma provides a wide range of bonded packing and configurations in the Novafix™ HPLC Cartridge System. This includes the most popular bonded packings on the market as well as packings for special applications. The Novafix Cartridge versatility of packings represents a great advantage over other cartridge systems that normally limit the range of packings to one or a few select packings.

- Packings of 3, 5 and 10 µm
- Lengths of 7.5, 15 and 25cm
- Different Internal diameters
- Packings of Tracer Excel, Tracer Extrasil, Lichrosorb, Lichrospher, Superspher.

## Guarantee

At Teknokroma, we guarantee the maximum quality of our products. This starts with quality in the mechanical components and finishes with final computerized quality control tests on each Novafix HPLC cartridge. Our quality controls ensure that you will receive only those cartridges which conform to the high quality demanded in our Novafix Cartridge specifications.

## Economy

To the criteria of maximum functionality and quality, we have also integrated the criterion of economy in the Novafix™ Cartridge System. The Novabond™ bonding and packing processes are rigidly controlled to produce superior yields of high-quality products. The Novabond™ processing makes NovafixNovafix™ Cartridges the most economical choice in the global HPLC market. This enables us to reduce the price even further with our "economy sets" of three NovafixNovafix™ Cartridges.

## Guard Cartridge

Function	Pkg	Cat.Nbr.
Silica	5 units	TR-015325
ODS	5 units	TR-015326
CN	5 units	TR-015327
NH <sub>2</sub>	5 units	TR-015328
SAX	5 units	TR-015329
SCX	5 units	TR-015330
C8	5 units	TR-015510
Diol	5 units	TR-015511
C <sub>6</sub> H <sub>5</sub>	5 units	TR-015512
C1	5 units	TR-015513
Carbohid.	5 units	TR-015331
Anions	5 units	TR-015335

## Kits and Accesories

Description	Cat.Nbr.
Holder: Connecting accessories : 2 units	TR-015323
Teflon seals: 10 units	TR-015324