

SEPARATION OF PESTICIDES

Column: **Meta.X5**, P/N TR-820232
Dimensions: 30m x 0.25mm x 0.25µm
Injection: 1 µL standard (1950 mg/L), split 1:20, 200°C
Carrier gas: He, 1mL/min
Oven program: 70°C(1min) @ 20°C/min to 150°C @ 3°C/min to 200°C
Detector: ECD, 250°C

Chromatogram provided by AINIA

<u>Peak Name</u>	<u>RT (min)</u>
1 Procolaz	18.99
2 Octacosane (Internal Standard)	21.02



TKG 1087