


OPTIMUN RESOLUTION IN SEMIVOLATILE COMPOUNDS ANALYSIS

Column: **Meta.X5**, 30m x 0.25mm x 0.5µm (P/N: TR-820532)
 Inj.: Splitless w/Surge: pulse 25psi @ 0.30min, 40ml/min @ 0.25min
 Inj. temp.: 250°C
 Oven temp.: 35°C (1min) to 280°C @ 25°C/min to 320°C (5min) @ 6°C/min
 Carrier gas: Helium, constant flow @ 1.2ml/min
 Det.: MS
 Transfer line temp.: 280°C
 Ionization mode: EI

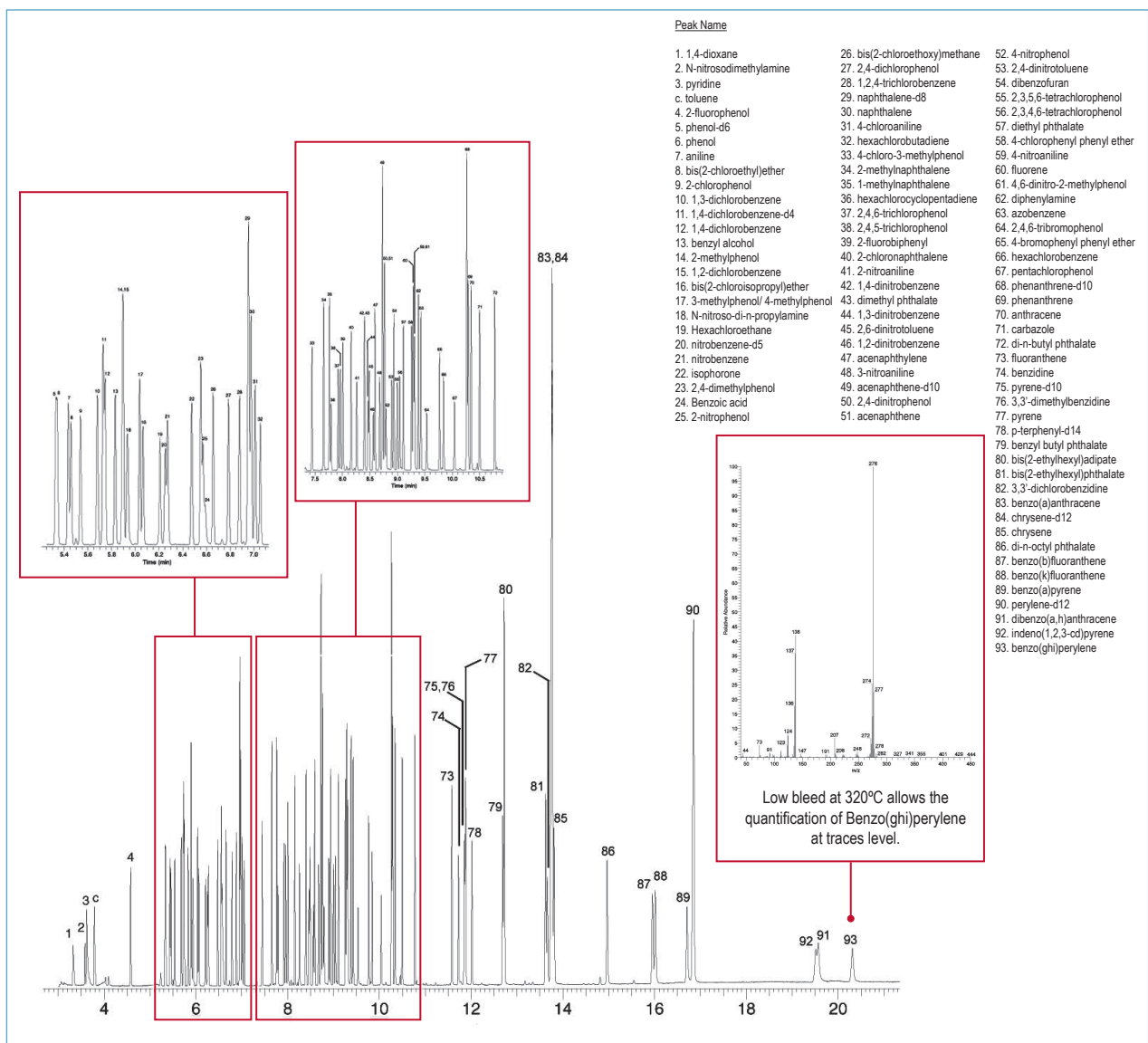

TKG 1258

www.teknokroma.com

Teknokroma S. Coop. C. Ltda. - Camí de Can Calders, 14 - 08173 Sant Cugat del Vallès (Barcelona) Spain
 Tel: +34) 936 748 800 Fax: +34) 936 752 405 e-mail: comercial@teknokroma.es